

3M Science.
Applied to Life.™

keydiagnostics

T: 02 8212 4074 F: 02 9423 6992
info@keydiagnostics.com.au
www.keydiagnostics.com.au
PO Box 1038, Gymea, NSW, 2227



The benefits are clear.

Get more than just results with 3M™ Allergen Testing.

**Simplify workflow and be confident in your food allergen testing process
with 3M™ Allergen Protein Rapid Kits and 3M™ Allergen Protein ELISA Kits.**

Testing for potentially harmful allergenic foods has become increasingly important to food and beverage processors and their customers. 3M™ Allergen Protein Test Kits provide qualitative and quantitative results and were designed for the quick and accurate detection of processed and unprocessed food allergen proteins.

Offering versatility across sample types, 3M Allergen Protein Test Kits can be used on environmental swabs, clean-in-place (CIP) final rinse water, raw ingredients and food samples. Now you can master the complex world of allergen testing. Streamline your process and get accurate results with 3M Allergen Protein Test Kits.

We've taken the hassle out of allergen testing. You see the results.



3M™ Allergen Protein Rapid Kits.



3M™ Allergen Protein ELISA Kits.

One tube process.

Sample type

Environmental swab
Clean-in-place (CIP) final rinse water
Raw ingredients
Finished food products



Add to 3M™ Extraction Buffer
Shake for 15 seconds



Pipette 100 µL to 3M™ LFD
Results in 11 minute ± 1 min.

Universal extraction protocol.


Sample type

(environmental swab, clean-in-place (CIP) final rinse water, raw ingredients, finished food products)

Perform one sample extraction



Test for multiple analytes

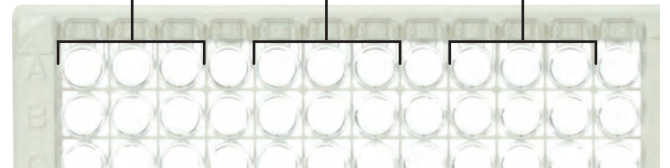

Walnut, pecan and crustacean proteins require a different extraction buffer, however the protocol is the same.

Complex testing made simple.

3M™ Environmental Collection Swabs and 3M™ Transfer Pipettes are available to complement 3M™ Allergen Protein Test Kits.



Ordering Information

Product Description	Catalog Number	Quantity
3M™ Environmental Collection Swab	ECS100	100 ea/pack
3M™ Transfer Pipette	TP100	100 ea/pack



3M Food Safety
3M Australia Pty Ltd
Bldg A, 1 Rivett Road
North Ryde NSW 2113

3M New Zealand Limited
94 Apollo Drive
Rosedale Auckland 0632

Phone 136 136
Web www.3M.com.au

Phone 0800 80 8182
Web www.3M.co.nz

See the complete portfolio.

3M.com.au/FoodSafety (Australia) or
3M.co.nz/FoodSafety (New Zealand)

3M is a trademarks of 3M.
Please recycle. © 3M 2019. All rights reserved.
AV011476031