

SpotCheck Plus™ | Food Safety Residue



SpotCheck Plus detects the presence of glucose and lactose on a surface as a rapid indicator of surface hygiene. Simply swab a surface, release the reagent, and if any residue containing glucose and/or lactose is present, the reagent will turn green. The more contamination present, the quicker the color change and darker the color. SpotCheck Plus instantly validates the hygiene of a surface, allowing immediate corrective action to be taken if necessary. If it turns green, it isn't clean!


Benefits:

- All-in-one device
- Easy-to-interpret color results
- No instrumentation necessary
- Results in 60 seconds
- Pre-moistened swab for reliable collection, recovery, and detection to ensure consistent results
- 100% recyclable
- 12 month shelf life at refrigerated temperatures (2-8°C)
- 4 week shelf life at room temperatures (21-25°C)

Sensitivity:

- Detects down to 2.5 μ moles D-glucose or 5.0 μ moles lactose in 60 seconds
- Disregard any color change after 2 minutes


Color Results:

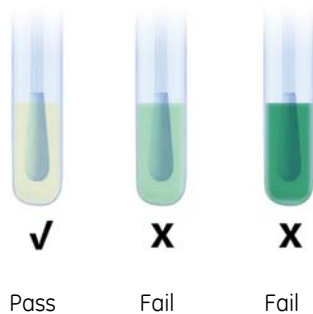

Find support documents, instructional videos, and more at www.hygiena.com

Catalog No.	Description	Quantity
HG-GL100	SpotCheck Plus	100

SpotCheck Plus Procedure:

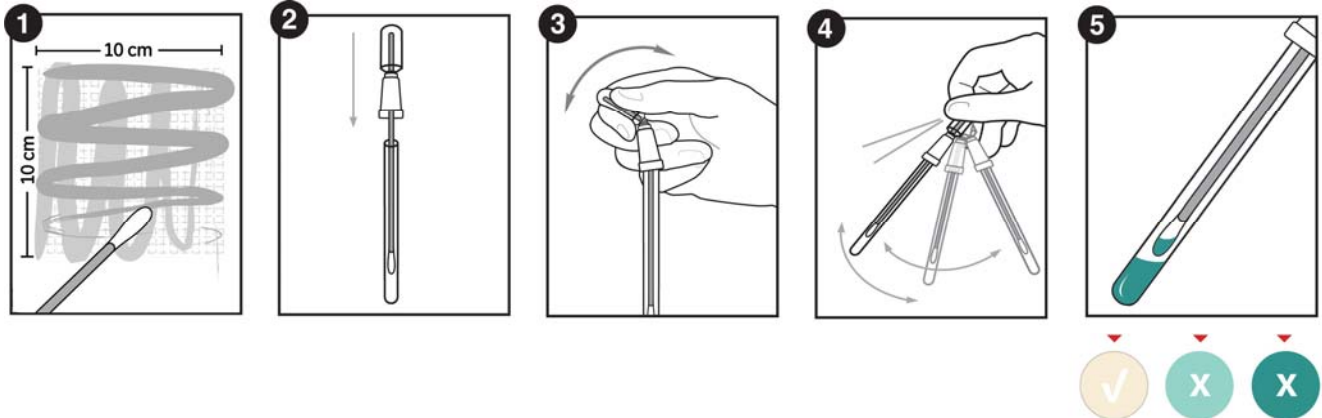

Related Products:



AllerSnap

AllerSnap is a colorimetric allergen prevention test for broad range allergen screening in 15-30 minutes. AllerSnap detects down to 3 µg protein.



ProClean

ProClean is a quick and easy-to-use device for monitoring the cleanliness of surfaces and safeguarding product quality. ProClean detects protein residues left on a surface after cleaning, giving results in 1-10 minutes without any required instrumentation.



SystemSURE Plus

SystemSURE Plus is used for ATP hygiene monitoring. This palm-sized instrument is easy to use, extremely sensitive and affordable, allowing users to quickly determine the cleaning efficiency and hygienic status of surfaces and water, ensure product quality, and reduce costs.

Catalog No.	Description	Quantity
HG-ALS-100	AllerSnap Allergen Prevention	100
HG-PRO100	ProClean Protein Residue Test	100
HG-SSPLUS	SystemSURE Plus ATP Monitoring System	1